



SMOBIO

Small Bio, Smart Tool

www.smobio.com

Product Information

ExcelDye™ 6X DNA Loading Dye, Tri-color

DL4000 5 ml × 2 (Tri-color, 6X)

Storage

4°C for 12 months

-20°C for 36 months

Working Reagent Preparation

1:6 dilution per DNA electrophoresis sample.

(1 μl loading dye for every 5 μl DNA sample)

For research use only
2019 ver. 2.1.1

P06-F11-A1

Description

SMOBIO's ExcelDye™ 6X DNA Loading dye (Tri-Color) is a pre-mixed buffer for tracking DNA samples during electrophoresis on agarose or polyacrylamide gels. It contains three dyes (Xylene cyanol FF, Bromophenol blue and Orange G) which help with the tracking of DNA migration. The Xylene cyanol FF, Bromophenol blue and Orange G migrate at approximately 800 bp, 150 bp and 30 bp on a standard 2% TAE agarose gel (4,000, 500 bp and 50 bp on 1% TAE agarose gel). The included glycerol helps keep the DNA at the bottom of the well, and the presence of EDTA chelates any divalent metal ions to prevent the process of a metal-dependent nuclease.

Composition

0.03% Xylene cyanol FF
0.03% Bromophenol blue
0.15% Orange G
10 mM Tris-HCl (pH 8.0)
60% glycerol
60 mM EDTA





SMOBIO
Small Bio, Smart Tool

FluoroBand DNA Ladders

