



**SMOBio**<sup>®</sup>

Small Bio, Smart Tool

[www.smobio.com](http://www.smobio.com)

## Product Information

ExcelTaq™ series

### Klen-Taq DNA Polymerase

---

#### TK1000 500 units

Klen-Taq DNA Polymerase (5 U/μl)	100 μl
10X Klen Buffer	1.2 ml

---

### Storage

-20°C for 24 months

### Features

- 5'→3' DNA polymerase activity
- 3'→5' exonuclease activity (proofreading)
- 4X fidelity as compared to *Taq* DNA polymerase
- Thermo-stable: up to 98°C during PCR denaturing step
- Robust PCR performance, resistant to variance in PCR conditions

## Description

The ExcelTaq™ Klen-Taq DNA Polymerase is a specially blended enzyme mix containing KlenTaq-1 DNA polymerase (a 5'-exo-minus, N-terminal deletion of *Taq* DNA polymerase) and a small amount of a proofreading DNA polymerase. This unique blending helps to improve the fidelity, yield and processivity of the resultant PCR process. Klen-Taq is also highly robust, showing high tolerance of varying concentrations of  $Mg^{2+}$ ; it is highly thermostable and has four times the fidelity compared to *Taq* DNA polymerase.

The ExcelTaq™ Klen-Taq DNA Polymerase is ideal for DNA amplifications 0.5-5 kb in length on genomic DNA, and up to 10 kb on less complex templates.

## **ExcelTaq™ Klen-Taq DNA Polymerase mixture**

Klen-Taq DNA Polymerase	5 U/μl
Proofreading DNA polymerase	Trace

### **Unit Definition**

One unit is defined as the amount of enzyme that will incorporate 10 nmol of dNTP into an acid-insoluble material in 30 minutes at 74°C.

### **Storage Buffer**

40 mM Tris-HCl (pH 7.5), 50 mM KCl, 0.1 mM EDTA, 25 mM (NH<sub>4</sub>)<sub>2</sub>SO<sub>4</sub>, 5.0 mM 2-mercaptoethanol, stabilizer, and 50% (v/v) glycerol

### **10X Klen Buffer**

400 mM Tricine-KOH (pH 9.2), 150 mM KOAc, 35 mM Mg (OAc)<sub>2</sub>, and 750 μg/ml BSA

## Recommended PCR Condition

Template	1 – 150 ng
Forward primer	0.1 – 0.5 $\mu$ M
Reverse primer	0.1 – 0.5 $\mu$ M
10X Klen Buffer	5 $\mu$ l
dNTPs	0.2 mM (each)
Klen-Taq enzyme	0.5 $\mu$ l (2.5 U)
ddH <sub>2</sub> O	to 50 $\mu$ l
<hr/>	
Total volume	50 $\mu$ l

## Recommended PCR Program

94°C	2 min	} 25 ~ 40 cycles
94°C	30 sec	
50~68°C*	30 sec	
68°C	30 sec/kb	
68°C	1 min	

\*Optimal PCR condition varies according to primers' thermodynamic properties.

## **Quality Control**

### **Functional Testing**

ExcelTaq™ Klen-Taq DNA Polymerase is tested for performance in the polymerase chain reaction (PCR) using 0.5 µl of enzyme to amplify a 665 bp target from 1 ng of tested plasmid DNA. The resulting PCR product is visualized as a single band on an ethidium bromide-stained agarose gel.

### **Nuclease Assay**

No contaminating endonuclease or exonuclease activity was detected using pUC19 incubated with ExcelTaq™ Klen-Taq DNA Polymerase for 4 hours at 37°C.

### **Residual Nucleotides Assay**

No contaminating residual nucleotides were detected from purified ExcelTaq™ Klen-Taq DNA Polymerase by PCR assay.

## **Other Information**

SMOBIO Technology, Inc. claims all warranties with respect to this document, expressed or implied, including but not limited to those of merchantability or fitness for a particular purpose. In no event shall SMOBIO Technology, Inc. be liable, whether in contract, tort, warranty, or under any statute or any other basis for special, incidental, indirect, punitive, multiple or consequential damages in connection with or arising from this document, including but not limited to the use thereof.

Caution: Not intended for human or animal diagnostic or therapeutic uses.

## Related Products

CK1000	Champion E. coli Transformation Kit
DM1100	ExcelBand 50 bp DNA Ladder, 500 $\mu$ l
DM2100	ExcelBand 100 bp DNA Ladder, 500 $\mu$ l
DM2300	ExcelBand 100 bp+3K DNA Ladder, 500 $\mu$ l
DM3100	ExcelBand 1 KB (0.25-10 kb) DNA Ladder, 500 $\mu$ l
NS1000	FluoroVue Nucleic Acid Gel Stain (10,000X), 500 $\mu$ l
RP1000	ExcelRT Reverse Transcriptase, 20,000 U
TF1000	SMO-HiFi DNA Polymerase, 100 U
TF3000	G-HiFi DNA Polymerase, 100 U
TP1000	ExcelTaq Taq DNA Polymerase, 500 U
TP1200	ExcelTaq 5 $\times$ PCR Master Dye Mix, 200 RXN
TP1260	ExcelTaq 5 $\times$ Fluorescent PCR Master Mix, 200 RXN
TP2100	ExcelTaq Blood Direct PCR Master Mix Kit, 200 RXN
TP5000	ExcelTaq Hot Start II DNA Polymerase, 500 U
TQ1110	ExcelTaq 2 $\times$ Q-PCR Master Mix (SYBR, ROX), 200 RXN
VE0100	B-BOX Blue Light LED epi-illuminator, AC 100-240V, 50/60Hz



## B-BOX™ Blue Light LED epi-illuminator

The latest version of the manual can be downloaded from [www.smobio.com/shop](http://www.smobio.com/shop).

For research use only  
2023 ver. 2.3.0